

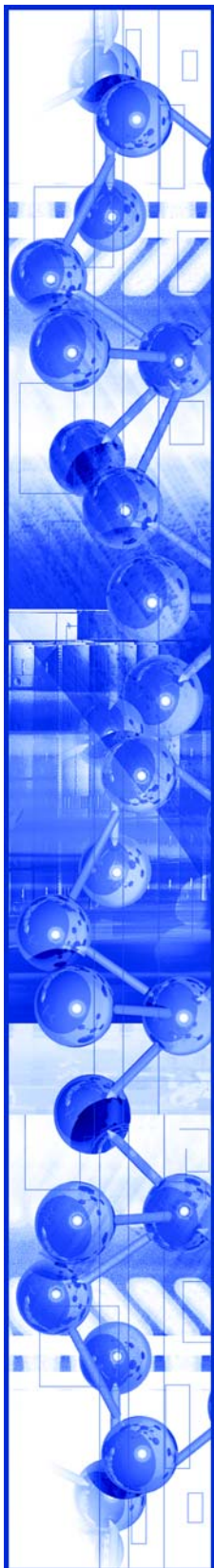
Rapid Western Blotting Kit

<u>Code</u>	<u>Description</u>	<u>Size</u>
N790-KIT	Rapid Western Blotting Kit-Mouse-SDT (<i>semi-dry transfer</i>)	1 Kit
N791-KIT	Rapid Western Blotting Kit-Rabbit-SDT (<i>semi-dry transfer</i>)	1 Kit
1B1322-KIT	Rapid Western Blotting Kit-Goat-SDT (<i>semi-dry transfer</i>)	1 Kit
1B1324-KIT	Rapid Western Blotting Kit-Rat-SDT (<i>semi-dry transfer</i>)	1 Kit
N792-KIT	Rapid Western Blotting Kit-Mouse-WT (<i>wet transfer</i>)	1 Kit
N793-KIT	Rapid Western Blotting Kit-Rabbit-WT (<i>wet transfer</i>)	1 Kit
1B1323-KIT	Rapid Western Blotting Kit-Goat-WT (<i>wet transfer</i>)	1 Kit
1B1325-KIT	Rapid Western Blotting Kit-Rat-WT (<i>wet transfer</i>)	1 Kit
	<i>Each kit includes</i> <i>Rapid Transfer Buffer, 10X, 150 ml</i> <i>RapidBlock™ Solution, 10X, 15 ml</i> <i>Rapid Antibody Diluent, 10X, 15 ml</i> <i>Rapid Wash Solution, 20X, 75 ml</i> <i>Rapid Western Anti-Mouse HRP, Rapid Western Anti-Rabbit HRP, Rapid Western Anti-Goat HRP, or Rapid Western Anti-Rat HRP, 30ul</i> <i>Sufficient material for 15 blots</i>	
N868-KIT	RAPID WESTERN BLOTTING KIT – SAMPLE <i>Kit includes</i> <i>Rapid Transfer Buffer, 10X, 2 x100 ml</i> <i>RapidBlock™ Solution, 10X, 2 ml</i> <i>Rapid Antibody Diluent, 10X, 2 ml</i> <i>Rapid Wash Solution, 20X, 6 ml</i> <i>Rapid Western Anti-Mouse HRP, 5 µl</i> <i>Rapid Western Anti-Rabbit HRP, 5 µl</i>	1 Kit

General Information:

AMRESKO's Rapid Western Blotting Kit streamlines each step of the Western Blotting procedure to reduce blotting time to about an hour. These versatile kits include all reagents except for user supplied primary antibodies and chemiluminescent detection reagents. No expensive transfer equipment is needed. The kit directly reduces the time for transfer and blocking steps and combines the primary and secondary incubations into a single step for additional time savings. Times and features for each step include:

- Transfer from gel to membrane with Rapid Transfer Buffer:
 - 10-20 minutes
 - Compatible with both wet and semi-dry apparatus
 - PVDF and nitrocellulose membranes
- Membrane blocking with RapidBlock™:
 - 5 minutes blocking time
 - Compatible with both PVDF and nitrocellulose membranes

**General Information:**

- Combined primary and secondary antibody incubation:
 - Complete in 30-45 minutes
 - Available with either anti-mouse or anti-rabbit HRP conjugated secondary antibodies

All reagents (excepting secondary antibodies) are protein-free formulations that minimize cross-reactivity and enhance antigen availability. Rapid Western Blotting Kit may be used with either PVDF or nitrocellulose membranes.

Storage/Stability:

All components of the Rapid Western Blotting Kit (**EXCEPT SECONDARY ANTIBODY**) are stable as concentrated or 1X solutions for 1 year at room temperature (18-25°C).

Rapid Western HRP-conjugated antibodies should be aliquoted and stored at -20°C for up to 1 year. Freeze-thaw cycles should be avoided. Antibodies are stable several weeks if kept undiluted at 2-8°C.

Application Disclaimer

*For research use only.
Not for therapeutic or diagnostic use.*



Protocol:

Reagents not included:

Deionized water

Chemiluminescent detection reagents

Notes:

- Rapid Western Blotting Kit is suitable for use with either PVDF or nitrocellulose membranes
- As with any immunoblotting procedure, antibody concentrations may need to be optimized empirically.
- Rapid Western Blot Secondary Antibody should be diluted immediately before use.

1. Transfer to Membrane

a. Semi-dry Transfer Apparatus

- Prepare a sufficient volume of 1X Rapid Transfer Buffer by diluting 1 part 10X Rapid Transfer Buffer into 9 parts distilled, deionized water.
- Gel preparation:
 - Wash gels in dH₂O for 5 minutes
 - Equilibrate in 1X Rapid Transfer Buffer for 5 minutes.
- Membrane preparation:
 - PVDF membranes: Prior to use, pre-wet in 100% methanol, rinse briefly in dH₂O, then equilibrate in 1X Rapid Transfer Buffer for a minimum of 5 minutes.
 - Nitrocellulose membranes: Equilibrate in transfer buffer prior to use.
- Assemble gel sandwich and semi-dry transfer unit according to manufacturer's instructions.
- Transfer at 25V for 10 minutes at room temperature.

b. Wet Transfer Apparatus

- Prepare a sufficient volume of 1X Rapid Transfer Buffer by diluting 1 part 10X Rapid Transfer Buffer into 9 parts distilled, deionized water.
- Membrane preparation:
 - PVDF membranes: Prior to use, pre-wet in 100% methanol, rinse briefly in dH₂O, then equilibrate in 1X Rapid Transfer Buffer for a minimum of 5 minutes.
 - Nitrocellulose membranes: Equilibrate in transfer buffer for a minimum of 5 minutes prior to use.

- Assemble gel sandwich and wet transfer unit according to manufacturer's instructions.
- Transfer at 75V for 20 minutes at room temperature.

2. Blocking

- Prepare a sufficient volume of 1X RapidBlock™ by diluting 1 part 10X RapidBlock™ Solution into 9 parts distilled, deionized water.
- Submerge membrane in 1X RapidBlock™ and incubate for 5 minutes at room temperature with agitation.

3. Antibody Incubation

- Prepare a sufficient volume (typically 5-10 ml per blot) of 1X Rapid Antibody Diluent by diluting 1 part 10X Rapid Antibody Diluent into 9 parts distilled, deionized water.
- Dilute primary antibody as recommended by the supplier into 1X Rapid Antibody Diluent.
- Dilute Rapid Western Secondary Antibody, HRP conjugated (N801-30UL, N802-30UL, 1B1311-30UL, or 1B1312-30UL into 1X Rapid Antibody Diluent solution containing primary antibody.

Example:

1X Rapid Blot Antibody Diluent	Primary Antibody	Rapid Blot Secondary Antibody, HRP Conjugated
10 ml	Supplier's recommendation	2ul

- Incubate the blot in antibody solution for 30-45 minutes at room temperature with agitation.

4. Washing

- Prepare a sufficient volume of 1X Rapid Wash Solution by diluting 1 part 10X Rapid Wash Solution into 9 parts distilled, deionized water.
- Wash the blot three times for 5 minutes each in 20ml 1X Rapid Wash Solution.

5. Detection

Visualize bands with enhanced chemiluminescent substrates for HRP linked reporter enzymes, such as AMRESKO's VisiGlo™ HRP Plus (N219-KIT).



Frequently Asked Questions

Questions

Answers

Why is the signal on my blot weak?

1. Primary antibody binds to the antigen with low affinity. Increase the concentration of primary antibody or the antibody incubation time.
2. Insufficient protein loaded on the gel. Load the SDS-PAGE gel with more protein.
3. Insufficient transfer of protein to the blotting membrane. Optimize transfer conditions and assess efficiency with Ponceau S or ProAct™ Membrane Stain.

Why is the background high on my blot?

1. Primary antibody concentration is too high. Reduce the amount of antibody used or the antibody incubation period.
2. Inadequate blocking of the membrane. Increase blocking time.
3. Methanol was used in the transfer buffer. Use only Rapid Transfer Buffer supplied with the kit or another methanol-free transfer buffer.
4. Insufficient washing. Increase the number and/or length of wash steps.

Why are there white bands on a black background on my blot?

1. Negative staining indicates antibody concentration is too high. Optimize antibody concentration and incubation time.

Can the Rapid Western Kits be used with phospho-specific antibodies?

Yes. The kits are compatible with phospho-specific antibodies.

Can I use a different transfer method in combination with the blotting kit?

If using a transfer buffer other than the Rapid Transfer Buffer, 10X supplied with the kit, you must use a methanol-free replacement. Methanol in the transfer buffer increases background on the blot.

Can I use Rapid Transfer Buffer with the iBlot® Dry Blotting System or other similar system?

Rapid Transfer Buffer is not compatible with devices, such as the iBlot® Dry Blotting System.

Related Products

M317	Fluorescent SPRINT NEXT GEL®, 10% Solution
M318	Fluorescent SPRINT NEXT GEL®, 12.5% Solution
M282	Proact™ Membrane Stain
K793	Ponceau S Stain
N218	VisiGlo™ HRP Chemiluminescent Substrate Kit
N219	VisiGlo Plus™ HRP Chemiluminescent Substrate Kit
N552	Gentle Review™ Stripping Buffer
N656	UnDo™ X-Ray Film Background Reducer

ZY0513 Rev. 1 05/2011 ©Copyright 2010 by AMRESKO LLC. All Rights Reserved. AMRESKO® is a registered trademark of AMRESKO, LLC.



Corporate Headquarters

AMRESKO LLC.
6681 Cochran Road
Solon, OH 44139 USA
Tel: (440) 349-1199
Toll free: 1-800-448-4442 (US & Canada)
Fax: (440) 349-2441
e-mail: info@amresco-inc.com

To Order (USA & Canada)

www.amresco-inc.com
Toll free: 1-800-448-4442
Toll free fax: 1-800-326-3733
Fax: (440) 349-3255
e-mail: dom-sales@amresco-inc.com

International Orders:

Tel: (440) 349-1199
Fax: (440) 349-2441
e-mail: int-sales@amresco-inc.com

Technical Support

Toll free: 1-800-610-2789 (USA & Canada)
Fax: (440) 349-0235
e-mail: techinquiry@amresco-inc.com